

# LABINCO L746 6 position Magnetic Stirrer 6 x 5 ltr.



## Technical data

Number of stirring positions	6 positions in two rows of three.
Stirring quantity max.	6 x 5 ltr
Motor rating input	6 x 35W
Motor rating output	6 x 15 W
Speed display	analog
Speed range	50 – 2500 rpm
Stirring bar length max.	60 mm
Heat output	-
Heating rate	-
Heating temperature range	-
Heat control	-
Heat control accuracy	-
Speed control	stepless
Adjustable safety circuit min.	-
Adjustable safety circuit max.	-
Connection for ext. temperature sensor	-
Control accuracy with sensor	-
Temperature constancy in medium	-
Set-up plate material	Stainles steel
Set-up plate dimensions	495 x 360 mm
Dimensions (W x D x H)	510 x 405 x 145 mm
Weight	12 kg
Permissible ambient temperature	5 - 40 °C
Permissible relative humidity	80 %
Protection class according to DIN EN 60529	IP 42
RS 232 interface	-
Product number	74600
Voltage	220 - 230 V
Frequency	50/60 Hz
Power input	220 W

# LABINCO L746 6 position Magnetic Stirrer 6 x 5 ltr.



## Description;

The "LABINCO MODEL L746 " is an instrument that provides stirring in beaker glasses, erlenmeyers or other suitable containers.

Stainless steel top plate.

Stirring speed is stepless controlled between 50-2500 r.p.m.

Robust housing and spillage protected control panel are provided, as well as integral lattice rod support at the rear.

Line cord and Teflon coated magnetic stirring bar are included with this model, delivered by ; Labinco bv, Breda, The Netherlands.

## Operating instruction;

After unpacking, place unit on a suitable surface. Put line cord into a suitable power-line (see specifications) and switch on " power".

The speed control works stepless, turn the knob clockwise until it turns.

Off course you have to introduce a magnetic stirring bar into the container.

**REMARK:** Containers made from easely magnetisable materials may be disturb stirring action.

## LABINCO BV

P.O. BOX 3259 4800DG BREDA THE NETHERLANDS

ZINKSTRAAT 64 4823AC BREDA THE NETHERLANDS

TEL;0031(0)76-5417582

FAX;0031(0)76-5420173

[E-mail;info@labinco-bv.com](mailto:info@labinco-bv.com)

<http://www.labinco-bv.com>