# LABINCO LD844 PTFE/Teflon coated 4 position Hotplate Magnetic Stirrer 4 x 5 ltr. 350°C

## Technical data

Number of stirring positions	4 positions in two rows of two.
Stirring quantity max.	4 x 5 ltr
Motor rating input	4 x 35W
Motor rating output	4 x 15 W
Speed display	analog
Speed range	50 - 1250rpm
Stirring bar length max.	60 mm
Heat output	4 x 300 W
Heating rate	7 K/min
Heating temperature range	room temp.- 350°C
Heat control	digital
Heat control accuracy	1 ±K
Speed control	stepless
Adjustable safety circuit min.	yes
Adjustable safety circuit max.	yes
Connection for ext. temperature sensor	on request
Control accuracy with sensor	1 ±K
Temperature constancy in medium	1 ±K
Set-up plate material	PTFE/Teflon coated
Set-up plate dimensions	Ø 120 mm
Dimensions (W x D x H)	345 x 385 x 185 mm
Weight	10 kg
Permissible ambient temperature	5 - 40 °C
Permissible relative humidity	80 %
Protection class according to DIN EN 60529	IP 42
RS 485 interface	On request
Product number	84402
Voltage	220 - 230 V
Frequency	50/60 Hz
Power input	1350 W

# LABINCO MODEL LD-814/816/844/846

## Description

The "LABINCO" Model LD-814/844/816/846 is a 4 or 6 position instrument that provides separate or simultaneous heating and stirring in beaker glasses, erlenmeyers or other suitable containers. Aluminum (silum) hotplates 120 mm diam. (Stainless steel on request) Temperature from ambient up to 350 degr.C.

Stirring speed is stepless controlled between 50-1250 r.p.m.

**Attention!! Do not set the temperature higher than 300 degrees**

## Operating instructions

After unpacking, place the model on a suitable surface. Put the line cord into a suitable power-line (see specifications) and switch on "power" on the backside. The speed control works stepless.

For the temperature setting, press the button " ✓ " and then set the desired temperature using the up or down arrow until the desired temperature on the screen and then again on " ✓ ".

The controller now regulated the temperature at the set value .

## LABINCO BV

P.O.BOX 3259	4800DG	BREDA	THE NETHERLANDS
ZINKSTRAAT 64	4823AC	BREDA	THE NETHERLANDS
TEL;0031(0)76-5417582			FAX;0031(0)76-5420173
<a href="mailto:info@labinco-bv.com">E-mail;info@labinco-bv.com</a>			<a href="http://www.labinco-bv.com">http://www.labinco-bv.com</a>

### Temperature setting:

Press once " ✓ " en now set the desired temperature using the up and down arrows. When the desired temperature is set once again to the 'set' button, the display shows the plate temperature and the temperature will go to the preset value.

### Time setting:

Press the right button at the top and set the desired time using the up and down arrows. When the desired time is set once again to the ' ✓ ' button, the display shows the plate temperature. In order to switch on the timer, 3 sec. The top right button pressed until it left a red light goes on with the heat signal and the timer is now on. To deactivate the timer hold the top right button again for 3 seconds. Pressed until the light goes out.

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