


LABINCO LD14 Large Magnetic Stirrer 60 ltr.

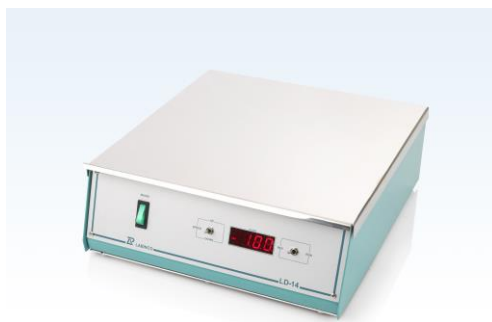

Technical data

Number of stirring positions	1
Stirring quantity max.	60 ltr
Motor rating input	50W
Motor rating output	20 W
Speed display	digital
Speed range	30 - 1000rpm
Stirring bar length max.	120 mm
Heat output	-
Heating rate	-
Heating temperature range	-
Heat control	-
Heat control accuracy	-
Speed control	stepless
Adjustable safety circuit min.	-
Adjustable safety circuit max.	-
Connection for ext. temperature sensor	-
Control accuracy with sensor	-
Temperature constancy in medium	-
Set-up plate material	Stainless steel
Set-up plate dimensions	320 x 320 mm
Dimensions (W x D x H)	320 x 340 x 105 mm
Weight	5.2 kg
Permissible ambient temperature	5 - 40 °C
Permissible relative humidity	80 %
Protection class according to DIN EN 60529	IP 42
RS 232 interface	-
Product number	14000
Voltage	100 - 230 V
Frequency	50/60 Hz
Power input	50 W

LABINCO LD14 Large Magnetic Stirrer 60 ltr.



Description

The Labinco model LD-14 is an instrument that provides stirring in beaker glasses, erlenmeyers or other suitable containers. *Stainless steel topplate 300 mm. *Stirring speed is stepless controlled between 30-999 RPM. *Robust housing and spillage protected control panel are provided, as well as integral lattice rod support at the rear. *Line cord and Teflon coated magnetic stirring bar are included with this model, delivered by Labinco bv, Breda, The Netherlands.

Operating instructions

After unpacking, place model LD-14 on a suitable surface. Put the line cord into a suitable power-line (see specifications). The speed control works stepless, push the speed-control knob up for speed till the stirrer starts rotating; the digital window will lit. The speed control can now be set for the required stirring action. Of course you have to introduce a magnetic stirring bar into the container.

Remark: Containers made from easily magnetizable materials may be disturb stirring action.

Labinco bv

P.O.BOX 3259	4800DG	BREDA	THE NETHERLANDS
ZINKSTRAAT 64	4823AC	BREDA	THE NETHERLANDS
TEL; 0031(0)76-5417582			FAX; 0031(0)76-5420173
E-mail;info@labinco-bv.com			http://www.labinco-bv.com