


# **NEW!! LABINCO LD208650 Rolling Bench 650 mm long rolls, with very strong motor up to 60 Kg, 600 Rpm and touchscreen!!**

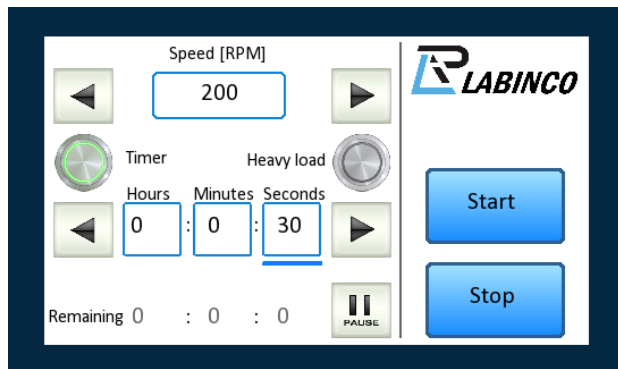


## **Technical data**

Number of rolling positions	1 up to 4
rolling quantity max.	60 kg
Motor rating input	250W
Motor rating output	175 W
Speed display	Digital
Speed range	1-600 rpm
Timer display	digital 1s. up to 999h 59m 59s
Heat output	-
Heating rate	-
Heating temperature range	-
Heat control	-
Heat control accuracy	-
Speed control	Digital
Adjustable safety circuit min.	-
Adjustable safety circuit max.	-
Connection for ext. temperature sensor	-
Control accuracy with sensor	-
Temperature constancy in medium	-
Set-up plate material	Steel with PE coated
Set-up dimensions	12 up to 2150 mm
Dimensions (W x D x H)	880 x 350 x 130 mm
Weight	11.5 kg
Permissible ambient temperature	5 - 50 °C
Permissible relative humidity	80 %
Protection class according to DIN EN 60529	IP 42
RS 232 interface	-
Product number	20865
Voltage	220 - 230 V
Frequency	50/60 Hz
Power input	300W

# **NEW!! LABINCO LD208 Rolling Bench with very strong motor up to 60 Kg, 600 Rpm and touchscreen!!**

## Description:

Rolling benches with standard 2 rolls  $\varnothing$  46MM, one roll is driven, extra rolls available on request.

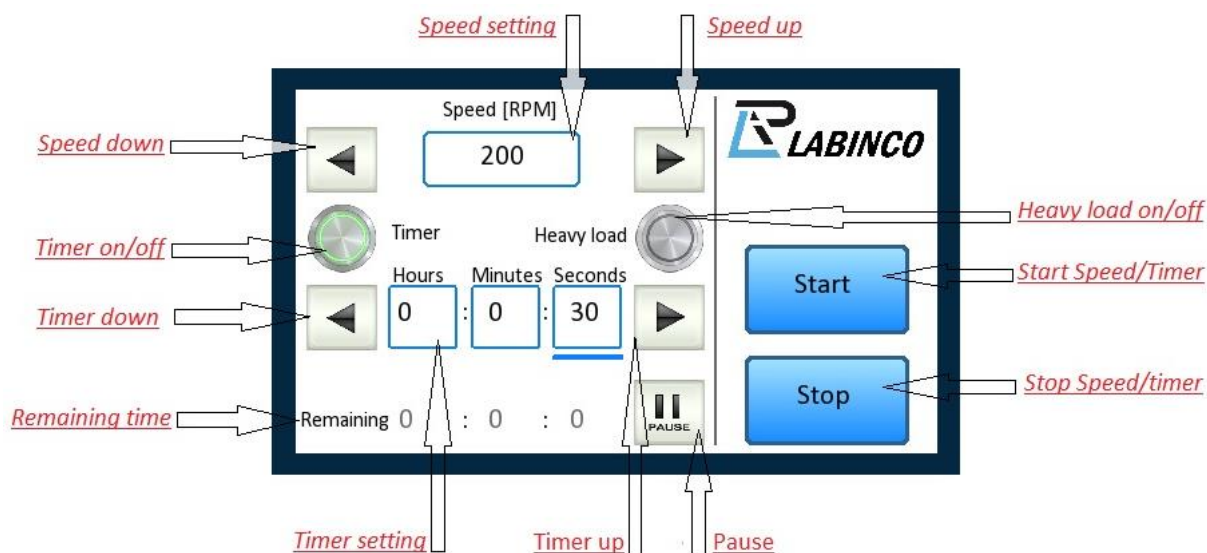

Operating instructions: Carefully unpack the unit and take off all packing materials. Choose a flat table or other suitable area for placing the Rolling bench. Check the model and catalog number indication at the rear. Check the line voltage in the room with the specified voltage on the label, this is standard 230 volts 50/60 HZ. Plug the line-plug into a suitable wall contact, with earth. Switch on the main switch. Pilot-lamp will light green if the bench is switched on.

This bench is equipped with standard 2 rolls diam.  $\varnothing$  46 mm, extra rolls available on request.

One roll is driven. You can easily place maximal 5 rolls on the rolling benches.

The first roll is driven, and the next roll is driven with the bottle or tube of the first roll, and the next also.... etc.

The minimal diam. is 12 mm up to ..... **But maximal 60 Kilogram**

Settings: You can easily set the speed or the time setting with the button "up" or "down" and if you are ready for start them, then push the start button, and if you will stop then, then push on the stop button. You can also for setting the speed or the timer function to touch the screen on the blue speed screen or timer screen, then the next screen will open for the setting.

## LABINCO BV

P.O.BOX 3259 4800DG BREDA THE NETHERLANDS

ZINKSTRAAT 64 4823AC BREDA THE NETHERLANDS

TEL;0031(0)76-5417582

FAX;0031(0)76-5420173

[E-mail; Info@labinco-bv.com](mailto:Info@labinco-bv.com)

<http://www.labinco-bv.com>